Exchangeable Thermoblocks for Thermomixers and Thermostats with IsoTherm-System

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eppendorf

Copyright© 2008 Eppendorf AG, Hamburg. No part of this publication may be reproduced without the prior permission of the copyright owner.

Trademarks

eppendorf is a registered trademark of Eppendorf AG, Hamburg, Germany.

Thermomixer is a trademark of Eppendorf AG, Hamburg, Germany.

Falcon is a registered trademark of Becton Dickinson, Franklin Lakes, USA.

Registered trademarks are not marked in all cases with TM or [®] in this manual

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1 User instructions

These instructions for use for exchangeable thermoblocks are a supplement to the operating manual for the thermomixer. They are not intended as a replacement for this operating manual.

Therefore, also read the operating manual of the corresponding unit before using the exchangeable thermoblock or the IsoTherm system for the first time. The current version of the operating manual can be found in the Internet at www.eppendorf.com.

2 Product description

2.1 Features

Exchangeable thermoblocks

The exchangeable thermoblock system for Thermomixer comfort/Thermomixer R and ThermoStat plus is used for efficient temperature control and mixing of all standard lab tubes.

The following exchangeable thermoblocks are available in combination with the IsoTherm System:

Exchangeable thermoblock	Tubes and plates	Use
Exchangeable thermoblock 0.5:	24 x 0.5 ml micro test tubes	<ul style="list-style-type: none">• Mixing and temperature control of the tubes.• Use Eppendorf IsoRack to reduce the heat exchange with the environment and maintain the temperature control.
Exchangeable thermoblock 1.5:	24 x 1.5 ml micro test tubes	
Exchangeable thermoblock 2.0:	24 x 2.0 ml micro test tubes	

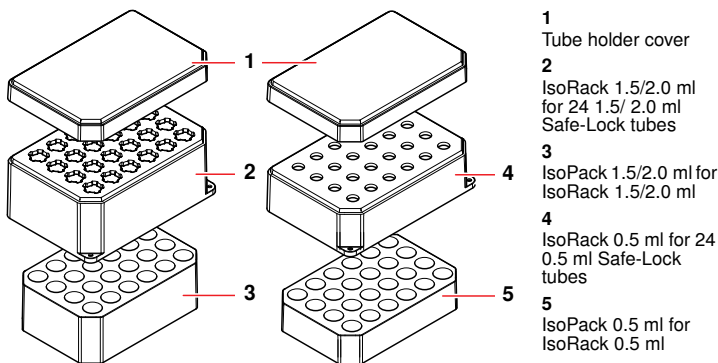
IsoTherm system

With the IsoTherm system samples can be cooled effectively and constantly over many hours and gently defrosted.

In addition, the micro test tube holder is used to assist in the filling of micro test tubes and in the provision, sorting, distribution, transport, storage and freezing of samples in micro test tubes.

The IsoTherm system includes the components IsoRack, IsoPack and IsoSafe.

IsoRack and IsoPack

Instructions for use

Component	Variants	Features
IsoRack (tube holder)	<ul style="list-style-type: none"> • for 0.5 ml Safe-Lock tubes, or • for 1.5/2.0 ml Safe-Lock tubes, each with cover. 	<ul style="list-style-type: none"> • base of an MTP • accepts 24 micro test tubes; numbered per row • temperature resistant between -200 and $+100$ °C • UV resistant and suitable for autoclaving • suitable for centrifuging in the MTP rotor with IsoRack adapter ¹ • stackable with or without cover • bore holes at the distances of the MTP grid, allowing for micro test tubes to be filled using multi-channel pipettes • cover fits twist-proof, locks to the tube holder • cover can also be used as base for cool pack and tube holder • floatable with cover (thawing in a water bath)
IsoPack (cool pack)	<ul style="list-style-type: none"> • for 0.5-ml IsoRack or • for 1.5/2.0-ml IsoRack 	both sizes are available in two versions each: <ul style="list-style-type: none"> • IsoPack white for 0 °C • IsoPack blue for -21 °C

¹ Generally the centrifuging stability depends on the centrifuge, rotor, rotor buckets, adapter, maximum rpm / maximum field of gravity (rcf), operating temperature, centrifugation times, density of the centrifuged liquid, organic solvents, accumulated total weight and proper operation.

3 Intended use

The exchangeable thermoblocks and the IsoTherm System are intended exclusively for indoor use. The exchangeable thermoblocks are used to hold tubes and plates, the IsoTherm system to cool tubes during sample preparation.

The exchangeable thermoblocks and the IsoTherm System must only be operated by appropriately qualified and trained specialist personnel. Carefully read the operating manual before starting work and familiarize yourself with the application.

4 Installation and operation

4.1 Mounting exchangeable thermoblocks

Warning! Injury or property damage from improperly attached exchangeable thermoblocks.

If the exchangeable thermoblocks are not attached properly, they can become detached from the device or break.

- ▶ Tighten the screws of the exchangeable thermoblocks fingertight.

Warning! Contamination due to opening tube cover.

Covers of micro test tubes can snap open involuntarily in the following cases, causing sample material to escape:

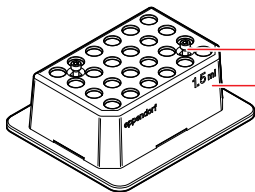
- high vapor pressure of the content
 - improperly sealed cover
 - damaged sealing lip
- ▶ Always check that micro test tubes have been sealed tightly before use.

Warning! Risk of burns from hot surfaces.

The exchangeable thermoblock can be very hot after heating and cause burns.

- ▶ Allow the heated exchangeable thermoblock to cool down completely before removing the exchangeable thermoblock.

1. Place the exchangeable thermoblock (2) on the thermosocket with the flat edge and the writing facing the front.
2. Tighten the captive screws (1).

4.2 Inserting tubes

- ▶ Press the tubes lightly into the holders of the thermoblocks.
To maximize temperature control, transfer the micro test tubes in IsoRacks to the thermoblock. The IsoRack reduces the heat exchange of the thermoblock with the environment. If necessary close the IsoRack cover.

4.3 Using the IsoTherm system

Caution! Damage of the IsoPack from sharp edges.

Sharp edged components can damage the film on the underside of the IsoPack.

- ▶ Do not use sharp edged components near the IsoPack.
-

Recommendations for precooling

- Cool pack 0 °C: approx. 2 hours (** freezer) or over night (* freezer)
- Cool pack -21 °C: approx. 4 hours at -35 °C (**** freezer, setting: rapid freeze)

Operation

1. Insert precooled cool box into the insulating box.
2. Place tube holder with samples onto the cool box.
3. Close insulating box with a cover.
This prevents the formation of frost.
The samples can be temperature controlled for several hours at 0 to -21 °C.

5 Maintenance

5.1 Cleaning

Clean the exchangeable thermoblocks and IsoTherm system components immediately if sample fluid enters the bore holes or the surface.

1. Clean the components with a mild soap solution. .
2. Rinse thoroughly with distilled water.

The IsoRack is UV resistant and can be autoclaved at 121 °C for 20 minutes. The tube holder and cover are resistant to alcohol.

Do not use organic solvents such as phenol, chloroform and acetone.

5.2 Disinfection / decontamination

Warning! Health hazard from contaminated accessories.

1. Follow the instructions in the decontamination certificate. It is available in PDF format on our homepage (www.eppendorf.com/decontamination).
 2. Decontaminate all the parts you want to dispatch.
 3. Include the fully completed decontamination certificate for returned goods in the package.
-

- ▶ Select a disinfection method complying with the statutory rules and regulations for your area of application. Use e.g. alcohol (ethanol, isopropanol) or alcohol-containing disinfectants.

6 Ordering information

Exchangeable thermoblocks

Order No. (International)	Order No. (North America)	Description
5361 000.015 5360 000.011 5362 000.019	022670506 022670522 022670549	Exchangeable thermoblock for microcentrifuge tubes complete with IsoTherm Rack and IsoTherm cool pack 0 °C 24 x 0.5 ml 24 x 1.5 ml 24 x 2.0 ml
5363 000.012	022670565	Exchangeable thermoblock for MTPs and deepwell plates with lid
5363 007.009	022670573	Adapter plate for 96 x 0.2 ml PCR tubes to insert in blocks for MTPs
5364 000.016	022670581	Exchangeable thermoblock for 24 micro test tubes 11 - 11.9 mm diam., height 30 - 76 mm
5366 000.013 5365 000.010	022670531 022670514	Exchangeable thermoblock for Falcon tubes for 8 x 15 ml for 4 x 50 ml
5367 000.017	022670557	Exchangeable thermoblock for 1.5 - 2 ml Cryo tubes
5368 000.010	022670590	Exchangeable thermoblock for slides holds 4 slides, for hybridization experiments
5368 000.100	022670786	Exchangeable thermoblock for slides DC holds 4 DualChip slides, with clamping plate
5352 100.007	022670603	Exchangeable thermoblock CombiBox for any test tubes interior 13.5 x 9.0 x 6.0 cm (W x D x H)
5355 000.011	022670107	Thermomixer comfort without exchangeable thermoblock

Instructions for use

Order No. (International)	Order No. (North America)	Description
5352 000.010	022670204	ThermoStat plus basic device without exchangeable thermoblock 100 - 240 V / 50 - 60 Hz

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Vaudaux-Eppendorf AG
Eppendorf AG · 22331 Hamburg · Germany · Tel: +49 40 538 01-0 · Fax: +49 40 538 01-556
Eppendorf North America, Inc. · 102 Motor Parkway, Suite 410 · Hauppauge, NY 11788-5178 · USA